

Hybridoma Cell Line

Quick Start Guide

Instructions

Preparations

Prepare concentrated liquid stocks by dissolving in cell culture grade water or PBS with gentle mixing or inversion. Sterile filter.

Storage

Liquid: stable up to 6 months at 4 to 8°C.

Powder: stable up to 12 months at -20°C.

Keep tightly sealed & protected from light.

Product	Test Concentrations	Notes & Recommendations
Zap-Hybridoma (100x Liquid) ¹	1x, 2x	DMEM/F12 is preferred classical medium
Cellastim	0.5, 1, 3 g/L	Broad effectiveness
Optiferrin	5, 10, 20 mg/L	Effective iron delivery
ITSE Animal-Free (100x Liquid)	1x, 2x	Reduces serum to 0.5% - 2%

1. Effective in Integra membrane partition systems.

For latest guidelines for use visit www.InVitria.com.

For technical support, please call your account manager or 1-800-916-8311. info@InVitria.com.

INVTRIA
Performance. Defined.